- DE Bedienungsanleitung
- GB Instruction Manual
- FR Mode d'emploi
- ES Manual de instrucciones

**BRESSER**  
Duolux

Art. No. 50-12000

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## RISK of physical injury!

Aids with sharp edges and tips are sometimes used with this device. Please store the device and all of its accessories and aids out of the reach of children. There is a risk of INJURY.

Children should only use this device under supervision. Keep packaging materials (plastic bags, rubber bands, etc.) away from children. There is a risk of SUFFOCATION.

The chemicals and liquids provided should be kept out of reach of children. Do not drink the chemicals! Hands should be washed thoroughly under running water after use. In case of accidental contact with the eyes or mouth rinse with water. Seek medical treatment for ailments arising from contact with the chemical substances and take the chemicals with you to the doctor.

This device contains electronic components that are powered by either a mains connection or batteries. Never leave a child unsupervised with this device. The device should only be used as per these instructions otherwise there is a serious RISK of ELECTRICAL SHOCK.

Batteries should be kept out of children's reach. When inserting batteries please ensure the polarity is correct. Leaking or damaged batteries can cause injury if they come into contact with the skin. If you need to handle such batteries please wear suitable safety gloves.

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Bresser GmbH

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## All parts (Fig. A):

- 1 16x WF eyepiece
- 2 5x WF eyepiece
- 3 Barlow lens
- 4 eyepiece barrel
- 5 microscope head
- 6 direct light adjustment (joystick)
- 7 LED lighting (direct light)
- 8 focussing wheel
- 9 combined colour filter and diaphragm wheel
- 10 battery compartment (base plate)
- 11 On/Off switch / dimmer
- 12 power connection
- 13 accessories tray
- 14 microscope table for transmitted light
- 15 microscope table for direct light
- 16 condenser disc for transmitted light
- 17 specimen plate for direct light
- 18 specimen holder clamps
- 19 objective
- 20 nosepiece
- 21 fixing screw
- 22 MicroOcular

## All parts (Fig. B):

- 23 MicroOcular with software
  - 24 Software initial screen
  - 25 MicroCut
  - 26 tweezers
  - 27 pipette
  - 28 preparation needles
  - 29 shrimp hatchery
  - 30 yeast
  - 31 Gum media
  - 32 sea salt
  - 33 shrimp eggs
  - 34 box of slides, covering glasses and permanent specimens
  - 35 mains plug
- 
- 36 cross table (optional)
  - 37 fastening screw
  - 38 longitudinal adjustment screw
  - 39 crosswise adjustment screw

## Index of contents

	Page
1. DuoLux Combi microscope	10
2. General/siting	10
3. Electric LED lighting with dimmer	10
4. Combined colour filter/diaphragm and wheel	10
5. Setting up the transmitted light table	10
6. Microscope settings (transmitted light)	10
7. Observation	11
8. Observation specimen - characteristics	11
9. Setting up the direct light table	11
10. Microscope settings (direct light)	11
11. Experiments (transmitted light)	11
12. Setup MicroOcular	12
13. Installing the software	12
14. Using the MicroOcular	13
15. Optional accessories	13
16. Care and maintenance	13
17. Disposal	13
18. Software troubleshooting	14
19. Technical data	14
20. EEC conformity explanation	14
Warranty & Service	25

## 1. DuoLux Combi Microscope

The DuoLux Combi Microscope is an entirely new product from Bresser. It combines the advantages of a direct and a transmitted light microscope with those of a handy combined microscope. The integrated changer slide in the microscope table allows you to convert the standard direct light table to transmitted light easily. This gives you an underlight workspace of about 38 mm diameter. The space between lens and table can be varied within a very wide range of around 45 mm so that high specimens can be examined, too. The direct lighting is a new development, too. It can be steered by joystick to the precise position on the specimen you require.

The lighting can be powered by either mains or battery. The combined colour filter and diaphragm wheel allows you to set optimal contrast in observing specimens using transmitted light. The integrated drawer for secure stowage of the accessories provided is another improvement.

## 2. General / siting

Before you begin assembling your microscope first select a suitable position for it. Make sure it stands on a stable base free of vibration.

Use of the electric lighting predicates mains power connection (220-230 V). (Use only the power connection/s supplied for safety reasons. Other such parts may not comply with the applicable technical specifications. We cannot accept any liability whatsoever for any damage due to the use of third-party power connection parts/plugs/sockets). Use the two alkali batteries type AA provided for battery operation. Insert them in the battery compartment in the base plate (Fig. A 10). Make sure the polarity is correct (shown on the battery compartment floor). Incorrect polarity will cause damage.

Accu batteries can be used but need to be charged using a special charger (neither is provided). The power supply unit provided is unsuitable for accu battery charging.

## 3. Electric LED lighting with dimmer

The microscope has two independently adjustable lighting units for

upper and lower lighting that can be controlled using the dimmer wheel (Fig. A 11). The transmitted light unit (Fig. A 9+14) is used for transparent specimens on holders/slides. To view solid non-transparent specimens use the direct light unit (Fig. A 7+15). Simultaneous use of both direct and transmitted light is only reasonable if the specimen is semi-transparent. This operating mode is not recommended for transmitted light specimens on slides as it may cause reflection on the slide.

## 4. Combined colour filter and diaphragm wheel

Use this (Fig. A 9), which is positioned in the slide tray under the microscope table, to adjust observation quality for transparent specimens. The finer details are better visible depending on colour and specimen. In direct light examination of e.g. transparent specimens coloured underlighting combined with white upper light may improve detail imaging.

Depending on the diaphragm opening used the appropriate light bundling can improve focus, focus depth, contrast and detail resolution.

## 5. Setting up the transmitted light table

Slide the changer to its stop in the table (as shown in Fig. A 14). Then insert the condenser disk (Fig. A 16) in the centre of table aperture.

## 6. Microscope settings (transmitted light)

The microscope view (Fig. A 1-5) is now adjusted for the first observation. First undo the screw (Fig. A 21) and turn the eyepiece barrel (Fig. A 4) until it reach a position comfortable for you.

You can now look through the lens, always beginning at the lowest magnification. Move the objective using the focussing wheel (Fig. A 8) to the top and turn the nosepiece (Fig. A 19) to the lowest magnification (4x). (The 4x objective is now positioned vertically downward).

### Note

Always move the objective right to the top before focussing. This prevents possible damage.

Insert the 5x eyepiece (Fig. A 2) in the Barlow lens (Fig. A 3). Make sure the Barlow lens is wholly inserted in the eyepiece barrel (Fig. A 4) and does not project.

## 7. Observation

Once you've set the microscope up with the right lighting and settings the principles below apply.

Start with simple observation at lowest magnification. Centring and adjustment of the specimen is thus made easier. The higher the magnification the more light is needed for good observation quality. Place a slide with a permanent specimen right under the lens on the microscope table and secure it in place with the clamps (Fig. A 18). Important. The specimen must be precisely sited in the centre of the transmitted lighting. Look through the eyepiece (Fig. A 1/2) and turn the focussing wheel (Fig. A 8) carefully until you get it right. Use the dimmer wheel (Fig. A 11) to set the brightness of the underlighting to view specimen details optimally. You can then set higher magnification by slowly pulling the Barlow lens (Fig. A 3) out of the eyepiece barrel (Fig. A 4). When nearly fully extracted magnification is nearly doubled.

For even higher magnification use the 16x eyepiece (Fig. A 1) and then turn the nosepiece (Fig. A 20) to a higher setting (10x/40x).

### Important note

Higher magnification need not necessarily lead to better results. This depends on the specimen.

### Please note

If magnification is changed (eyepiece or objective change, Barlow lens extraction) the focus must be re-adjusted using the focussing wheel (Fig. A 8). Be very cautious when doing this. If you turn the lens down too fast you may damage it and/or the slide.

## 8. Observation specimen - characteristics and preparation

### 8.1 Condition

Both transparent and non-transparent specimens can be examined with this microscope, which is a direct as well as transmitted light model. If opaque specimens are examined - such as small animals, plant parts, tissue, stone and so on - the light is reflected from the specimen through the objective and eyepiece, where it is magnified, to the eye (reflected light principle). If opaque specimens are examined the light from below goes through the specimen, objective and eyepiece to the eye and is magnified en route (direct light principle). Many small organisms of the water, plant parts and finest animal components have now from nature these transparent characteristic, other ones must be accordingly prepared. Is it that we make it by means of a pre-treatment or penetration with suitable materials (media) transparent or thus that we cut finest wafers off of them (hand cut, MicroCut) and these then examine. With these methods will us the following part make familiar.

### 8.2 Producing a thin specimen slide

As already stated, specimens for microscopic observation should always be sliced as thin as possible. A little wax or paraffin is needed to achieve the best results. A candle can be used for the purpose. The wax is put in a bowl and heated over a flame. The specimen is then dipped several times in the liquid wax. The wax is finally allowed to

harden. Use a MicroCut (Fig. B 25) or knife/scalpel (careful!!!) to slice the wax-coated specimen as thinly as possible. The slices are laid on slides and covered with a covering glass.

### 8.3 Making your own specimens

Place the specimen on the slide and add a drop of distilled water with a pipette (Fig. B 27) to the specimen (Fig. B 1).

Set a covering glass vertically at the edge of the drop so that the water runs along the glass edges (Fig. B 11). Then slowly place the covering slide atop the drop.

**Note: The gum media provided (Fig. B 31) is used in making permanent slides. Add it instead of distilled water. The medium hardens and the specimen is then permanently affixed to the slide.**

## 9. Setting up the direct light table

Pull the sliding change tray out of the table to its stop (as in Fig. A 15). Then insert the specimen plate for direct light (Fig. A 17) in the central table aperture.

## 10. Microscope settings (direct light)

The wide adjustment range of the microscope allows you to examine specimens up to a height of about 40 mm. The variable direct lighting can be precisely adjusted using the joystick (Fig. A 6) to the specimen. In examining transparent specimens the underlighting with its approximate diameter of 32 mm is an ideal aid to getting a really good view of the specimen. You can adjust both upper and lower lighting individually or together to optimally illuminate any specimen.

The best results in the direct light mode are achieved by combining the 5x eyepiece and the 4x objective. Any other combination increases magnification but reduces the field of visibility.

## 11. Experiments (transmitted light)

Once you're familiar with the microscope you can try the following experiments and view the results.

### 11.1 Newspaper print

#### Objects:

1. A small piece of paper from a newspaper with parts of a picture and some letters,
  2. a similar piece of paper from an illustrated magazine.
- Use your microscope at the lowest magnification and use the preparation of the daily paper. The letters seen are broken out, because the newspaper is printed on raw, inferior paper. Letters of the magazines appear smoother and more complete. The picture of the daily paper consists of many small points, which appear somewhat dirty. The pixels (raster points) of the magazine appear sharply.

### 11.2 Textile fibers

#### Objects and accessories:

1. Threads of different textiles: Cotton, line, wool, silk, Celanese, nylon etc.,
2. two needles.

Each thread is put on a glass slide and frayed with the help of the two needles. The threads are dampened and covered with a cover glass. The microscope is adjusted to a low magnification. Cotton fibres are of vegetable origin and look under the microscope like a flat, turned volume. The fibres are thicker and rounder at the edges than in the centre. Cotton fibres consist primary of long, collapsed tubes. Linen fibres are also of vegetable origin; they are round and run in straight

lines direction. The fibres shine like silk and exhibit countless swelling at the fibre pipe. Silk is of animal origin and consists of solid fibres of smaller diameter contrary to the hollow vegetable fibres. Each fibre is smooth and even moderate and has the appearance of a small glass rod. Wool fibres are also of animal origin; the surface consists of overlapping cases, which appear broken and wavy. If it is possible, compare wool fibres of different weaving mills. Consider thereby the different appearance of the fibres. Experts can determine from it the country of origin of wool. Celanese is like already the name says, artificially manufactured by a long chemical process. All fibres show hard, dark lines on the smooth, shining surface. The fibres crinkle after drying in the same condition. Observe the thing in common and differences.

### 11.3 How does bread mould develop?

**Object:** An old piece of bread.

The spores of the kind of mould, which strike our bread, are to be found everywhere in the atmosphere. Put bread on a slide and squirt carefully some water on it. Moisten bread only, don't wet it. Put the whole into a container with a screw-type cap and place it into a cabinet, into which only little light breaks in and which prevails it in a warm temperature. Within a short time the black bread mould forms. Regard the bread each day. At the first the mould shows up a white, shining consistence. Take it on a slide to observe it. The material turns out as a complicated thread mass, which forms the fungus body in its whole. One calls the whole mycelium. Each thread is a hypha. Soon some rhizoids arise, which embody the mould fungus with bread, in order thereby to water and nutrients for the growth of the mycelium to receive. In the course of time the rhizoid colours itself brownish. Vertically over this group hyphae grow like long slim stacks, which end in a tiny small, white ball. One calls the stack sporangiophores (carrier of the sporcap), the ball is a sporangium or a sporcap. Soon these balls accept a black color. Inside spores present mature. If now the sporcap breaks open, then it sets the spores free, which step now to air and infect other bread. With the naked eye you can recognize the mature sporcaps as tiny black marks. They are scattered on the mould fungus surface and give thus to the kind of fungus its name. There are however still different kinds of mould fungi. They can be pink, red, blue or green. Manufacture yourselves preparations of all stages of the bread mould.

### 11.4 Salt water shrimps

**Accessories:**

1. Yeast (Fig. B 30),
2. Gum media (Fig. B 31),
3. Sea salt (Fig. B 32),
4. Shrimp eggs (Fig. B 33),
5. Shrimp egg hatching plant (Fig. B 29).

#### 11.4.1 The lifecycle of the saltwater shrimp

The saltwater shrimp or *Artimia salina* to scientists has an unusual and interesting lifecycle. The female's eggs are hatched without any male shrimp having to fertilise them. The resultant baby shrimps are all female. Under unusual conditions such as when a swamp is drained the eggs may produce male shrimps. These males fertilise the female's eggs, resulting in a specific type of eggs. These are called winter eggs and have a thick shell as protection. They're pretty rugged and can survive the swamp or lake drying out causing the death of the entire shrimp population for up to a decade in a form of hibernation. The eggs hatch once the right ambient conditions again obtain. The eggs supplied (Fig B 33) are of this type.

#### 11.4.2 Hatching of the salt water shrimp

To hatch the shrimp it is essential to first have a saline solution suited to the shrimp's needs. Fill half a litre of rain- or freshwater in a container. Let it stand for about thirty hours. As water evaporates over time it's a good idea to have a second container of such water left standing for thirty-six hours. Once it's stood for this length of time pour half of the sea salt supplied into one of the containers and stir until it has dissolved. Then pour some of it into the shrimp's breeding plant (Fig. B 29). Add a few eggs and close the lid. Put it somewhere with plenty of light but not in the direct sun. The temperature should be approximately 25° C. The shrimps will hatch in two or three days at this temperature. Should any water evaporate during this time replace it from the second container.

#### 11.4.3 The saltwater shrimps under the microscope

What comes out of the egg is known as a nauplius larva. Use the pipette (Fig. B 27) to put some of them on a slide for examination. They will move in the solution using their hair like limbs. Remove a few daily from the container for examination under the microscope. If you do so and save the pictures made with the MicroOcular you will then have a seamless record of the shrimp's lifecycle. You can remove the upper lid of the shrimp's breeding plant and put the whole thing under the microscope. The larvae will mature in six to ten weeks depending on ambient temperature. You will soon have bred an entire generation of saltwater shrimps that constantly reproduce.

#### 11.4.4 Feeding your saltwater shrimps

To keep them alive saltwater shrimps must be fed occasionally. This must be done carefully as overfeeding causes the water to stagnate and poison the shrimps. Feeding is best done with dry powdered yeast (Fig. B 30). Give them a little every other day. If the water darkens this signifies it is stagnating. If so remove the shrimps and put them in a fresh saline solution.

**Note**  
Eggs and shrimps are unfit for human consumption!

**Note:** The MicroOcular only works without the Barlow lens supplied! The magnification setting is changed using the MicroOcular and has to be reset with the focussing wheel. To ensure proper lighting of the specimen using the MicroOcular working with batteries is not recommended. Please use the mains power adaptor supplied.

## 12. Setup MicroOcular

Remove the Barlow lens (Fig. A 3) with the currently used eyepiece from the tube (Fig. A 4) and put the MicroOcular (Fig. A 22) in place instead (Fig. A).

Please do not yet connect the MicroOcular to your PC. Please go through the following points step by step.

## 13. Installation and use of the software

### 13.1 Installation of the image processing software Photomizer SE

- 1 Insert the CD-ROM (Fig. B 23) supplied in your CD or DVD drive. The installation menu opens automatically. If it doesn't, please switch to Windows Explorer and select your CD drive. Start the file "setup.exe" by double-clicking it.
2. A choice of languages is shown. Choose your preference. Confirm it by clicking „OK“. At the „Welcome“ screen, click „Next>“.
4. In the next screen you will be asked for the „Destination Folder“. Confirm your input with „Next>“.

- The setup status screen shows. This advises current operation status with a progress bar. This may take a few minutes.
- The „Completed the Photomizer Setup Wizard“ screen shows. Click here on „Finish“.

#### Note

To use the MicroOcular long-term we recommend it always be connected to the same USB port.

### 13.2 MikrOkular driver installation

During Photomizer SE software installation the correct driver for your operating system will be installed automatically. No manual input is needed.

In rare cases the device may not be recognised by your computer. As a rule you need then only install the driver manually from the CD. If this fails please refer to the troubleshooting chapter that follows.

### 13.3 Troubleshooting: the device isn't recognised by the computer (USB)

- This occurs in rare cases whith some OEM Windows Vista or XP versions. The device is not defective. Updating the notebook or PC USB driver usually suffices.
- If not, proceed as below. (The maker cannot accept any liability whatsoever in these cases<sup>1</sup>).

Delete ghost devices. These are devices not actually currently connected to your computer. The reason for this is that Windows generates a system settings entry for each USB device ( e.g USB Stick) on each port. It then seeks that/those device/s each time it's started. This slows down USB device recognition and Windows booting and causes non-recognition of USB ports.

Cleaning up the system environment and deleting devices that are no longer in use helps here. To do so right click WORKPLACE (COMPUTER in Vista) and open the PROPERTIES menu. Then click the ADVANCED tab (ADVANCED SYSTEM SETTINGS in Vista) and then ENVIRONMENT VARIABLES. Choose the NEW option in ENVIRONMENT VARIABLES. Enter the following under VARIABLE NAME.

devmgr\_show\_nonpresent\_devices

Under VARIABLE VALUE enter „1“.

Confirm your input with OK and then reboot the computer.

Change to the device manager. In the VIEW menu activate the SHOW DEVICES OPTION. The ghost devices previously not shown will now be displayed in light grey. Check the various categories including USB, memory size, etc. Only delete entries for devices you no longer use.<sup>1</sup>

<sup>1</sup>Source: Microsoft Knowledge Base

<http://support.microsoft.com/kb/315539/en>

\*Not included

\*\*System requirements: Intel X86 processor with at least 2.0 GHz speed, 200 MB hard disc space, 512 MB RAM memory, Windows XP or Vista (32-bit) and a free USB port (not via a hub if possible).

## 14. Using the MicroOcular

### 14.1 Preparation

- Slide a specimen under your microscope and focus on it.
- Remove the eyepiece and Barlow lens from the eyepiece barrel and the dust cap from your MicroOcular and install same in the eyepiece barrel instead of the Barlow lens.
- Start your PC if you haven't yet and connect your MicroOcular to the USB port of your computer.

### 14.2 Showing and saving MikrOkular images on your PC

- Start the Photomizer SE Software (Fig. B 24).
- Click „Open camera“
- If you have more than one device connected up select the one you want from the choice offered. Click „SoC PC-Camera“ or „MikrOkular“. Omit this step if you only have a single device connected up.
- The camera image should now be visible on your screen. Focus the microscope image.
- Click „Capture“ to record an image. It will then be shown on the right.
- Click on it to select it and then click „Transfer image“.
- This takes you to the Photomizer SE software.
- File - Save as

### 14.3 The Photomizer SE software

If you need help in the „Photomizer SE“ programme click „?“ and then „Open help“. If you need further assistance please visit the maker's home page at [www.photomizer.net](http://www.photomizer.net)

## 15. Optional accessories

A cross table is available for the Bresser DuoLux as an optional accessory (Fig. B 36). Remove the specimen holder clamps to install it (Fig. A 18).

Use the screw (Fig. B 37) to fasten the cross table to the microscope table. Use the adjusting screws (Fig. B 38+39) to precisely position the specimen longitudinally (Fig. B 38) and crosswise (Fig. B 39).

## 16. Care and maintenance

Remove the device from it's energy source before cleaning (remove plug from socket / remove batteries). Lenses (eyepieces and/or object lenses) should be cleaned with a soft, lint-free cloth (e.g. microfibre) only. Do not use excessive pressure - this may scratch the lens. Clean the exterior of device with a dry cloth. Do not use cleaning fluids so as to avoid causing damage to electronic components. Protect the device from dust and moisture. Store the device in the bag supplied or in its original packaging. Batteries should be removed from the device if it is not going to be used for a long period of time.

## 17. Disposal

Dispose of the packaging material/s as legally required. Consult the local authority on the matter if necessary.

Do not dispose of electrical equipment in your ordinary refuse. The European guideline 2002/96/EU on Electronic and Electrical Equipment Waste and relevant laws applying to it require such used equipment to be separately collected and recycled in an environment-friendly manner. Empty batteries and accumulators must be disposed of separately. Information on disposing of all such equipment made after 01 June 2006 can be obtained from your local authority.

## 18. Software troubleshooting

Fault	Solution
No image visible	<ul style="list-style-type: none"><li>• Turn the lighting on</li><li>• Insert condenser lens (Fig. A 16) when viewing with the naked eye</li><li>• Re-focus</li></ul>
Image flickers (when viewing with the MicrOcular)	<ul style="list-style-type: none"><li>• reduce graphics card resolution</li></ul>
Software installation causes "not XP approved" to show	<ul style="list-style-type: none"><li>• confirm with &lt;OK&gt;</li></ul>

## 19. Technical data

MicrOcular system requirements: Intel Pentium 166 MHz or comparable or better, Windows 2000/ME/XP/Vista operating system, Min. 64 MB RAM, Min. 100 MB hard disc space

### Magnification table

Eyepiece	Objective	Magnification	with Barlow lens
5x	4x	20x	40x
5x	10x	50x	100x
5x	40x	200x	400x
16x	4x	64x	128x
16x	10x	160x	320x
16x	40x	640x	1280x

### Special aspects

Change table for direct and transmitted lighting. Extremely long adjustment range (45 mm) for viewing specimens up to max. 40 mm in direct lighting. Direct light table with transparent work space of about 32 mm diameter. Economical LED lighting for direct and transmitted lighting, combined or separate. Brightness for direct and transmitted lighting separately adjustable. Direct lighting adjustable by joystick. Combined colour filter and diaphragm wheel for both direct and transmitted lighting modes. Inbuilt tray for accessories. Battery or mains powered lighting.

## 20. EEC conformity explanation

Bresser GmbH, resident in 46414 Rhede/Westf., Gutenbergstr. 2, Germany, explains the agreement with in the following specified EEC guidelines for this product:

EN 55011/A2:2007  
EN 61000-6-1:2007  
EN 61000-3-2:2006  
EN 61000-3-3/A2:2005

Product description: BRESSER DuoLux  
Model: Biological Monocular Microscope

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## GB WARRANTY & SERVICE

The regular guarantee period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required. You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at [www.bresser.de/warranty\\_terms](http://www.bresser.de/warranty_terms).

Would you like detailed instructions for this product in a particular language? Then visit our website via the link below (QR code) for available versions.

Alternatively you can also send an email to [manuals@bresser.de](mailto:manuals@bresser.de) or leave a message on +49 (0) 28 72 - 80 74-220\*. Please always state your name, precise address, a valid phone number and email address, as well as the article number and name.

\*Number charged at local rates in Germany (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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