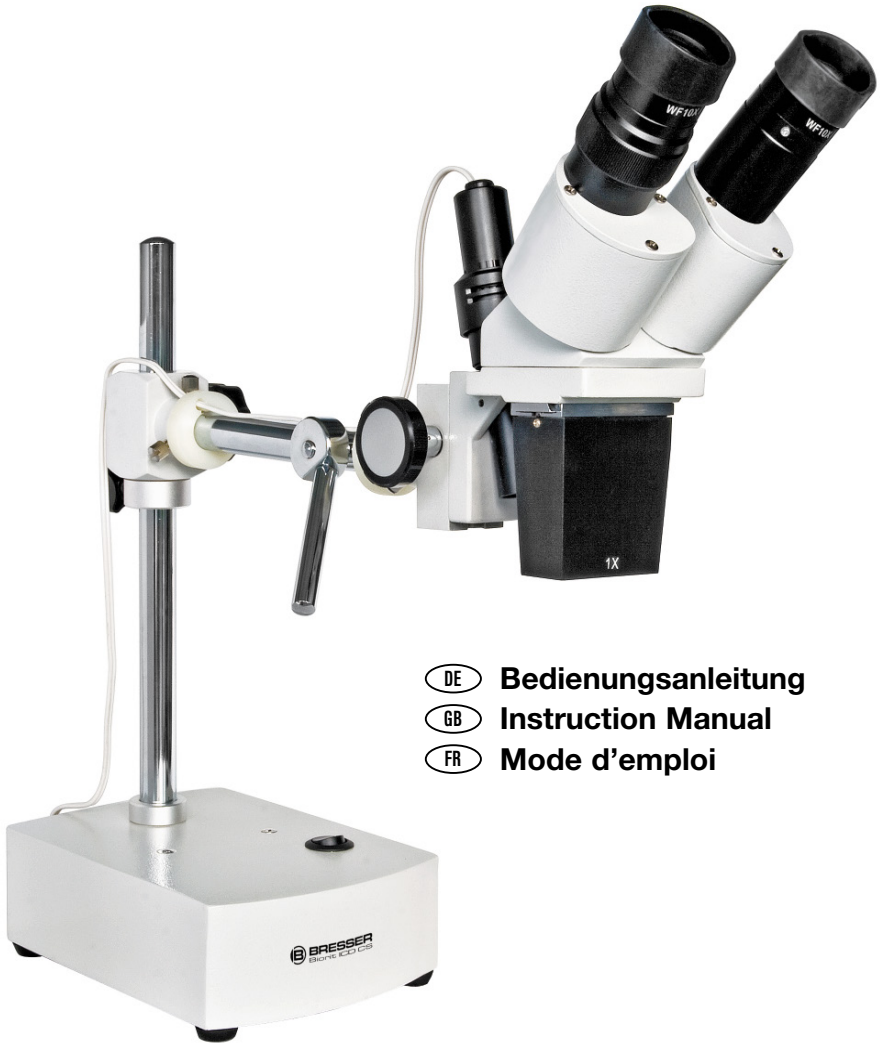
BRESSER®

- DE** Bedienungsanleitung
- GB** Instruction Manual
- FR** Mode d'emploi

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Telefon:.....Kaufdatum:

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CAUTION!

To work with this microscope, sharp and pointed aids are being used. Please take care that this microscope and its accessories are stored at a place out of reach of children. Let children only work with this microscope under an adult's supervision!

Keep packing material (plastic bags etc.) away from children!

1. Parts of the microscope

- | | |
|---------------------------|---------------------|
| ① Eyepieces | ⑧ Panhandle |
| ② Diopter adjustment ring | ⑨ Focusing wheel |
| ③ Illuminator | ⑩ Objective |
| ④ cable holder ring (a/b) | A Eyepieces, |
| ⑤ Vertical microscope arm | B 2 Rubber eyecups, |
| ⑥ Base | C Replacement lamp |
| ⑦ Power switch | |

Suitable location for observation:

Dry and dust free, room temperature between -5 and +40°C.

Position your device so that it can be disconnected from the power supply at any time. The wall socket should always be located near the device and be easily accessible, since the plug on the power cord serves as a disconnecting device for the power supply.

2. Assembly

- I. Place the base (6) on a stable surface.
- II. Place the first cable holder ring on the horizontal microscope arm.
- III. Place the second cable holder ring on the vertical microscope head pole.
- IV. Fix both arms together, using the handle (8) with the spring lock washer.
- V. Push the illuminator (3) through the hole opening in the microscope head.

3. Observation

Turn on the microscope using the power switch (7).

Now, look through the eyepieces (1) of the binocular viewing and change the spacing between the eyepieces so that you get a round, clear field of view. The definition is adjusted by slowly turning the focusing wheel (9).

If you have to observe a very large or very small object, you can adjust the microscope. Loosen the holding screw on the side and move appropriately the microscope head on its guidance. If you have found the correct position (distance of 22 cm from the object), tighten the holding screw again.

4. Diopter adjustment

Turn the left diopter adjustment ring (2) on the left eyepiece to the central position (silver marking). Shut your left eye and look with the right eye through the right eyepiece. Now focus with the focusing wheel (9), focus correctly for your right eye. Now, look with your left eye through the left eyepiece. If the image isn't sharp, adjust the focus for your left eye by turning the diopter adjustment ring. Now, the microscope is correctly adjusted.

5. Exchange the Lamp

Make sure that the microscope is separated from power outlet. Let the lamp cool down, if necessary, since lamps can become very hot in operation.

- I. Please unscrew the lamp housing set screw.
- II. Pull off the lamp housing.
- III. Now exchange the bulb.
- IV. Replace the lamp housing and fix it with the set screw.

6. Care and maintenance

Cleaning the lenses: Sweep the lens with a lens tissue or soft fabric immersed with a small amount of alcohol (e.g. isopropanol).

Cleaning the painted parts: The dust on the painted parts can be removed by cleaning cloth. For the grease spots, the gauze immersed slightly with aviation gasoline is recommended. Do not use organic solvents such as alcohol, ether or other thinner etc. for cleaning the painted parts or plastic components. Your microscope is a high-quality optical instrument. Avoid disassembling the microscope. Improper disassembling affects the high utility of the microscope. Moreover, the warranty expires. Cover the microscope with the dust protective cover and place it there, where it is dry and dust-free. We suggest the storage of all objectives and eyepieces in a closed container with drying agent.

7. Technical data

Binocular viewing

Eyepiece: WF 10x for 10x magnification, WF 20x (not included) for 20x magnification

Objective: 1x

Illumination: 12 V, 10 W (bulb)

Fuse: 0.5 A, 250 V

Accessories: Spare lamp, 2 rubber eyecups, AC adapter, dust protecting cover, instruction manual.

8. Only for EU countries

Do not dispose of electric equipment together with household waste material! In observance of European Directive 2002/96/EC on waste electrical and electronic equipment (WEEE) and its implementation in accordance with national law, electric equipment that have reached the end of its life must be collected separately and returned to an environmentally compatible recycling facility. Discharged batteries and damaged re-chargeable batteries must be disposed of at special battery collection points. Information is available from your local disposal agent or local authority regarding the disposal of devices or batteries.

9. Declaration of Conformity

Bresser GmbH has formed a „Conformity Declaration“ in line with the applicable guidelines. This declaration is available for review.

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10. Warranty

The period of warranty is 2 years, beginning on the day of purchase. Please keep the cash receipt as evidence of purchase. Devices which become defective during the warranty period can be returned to the dealer where the device was bought. The repaired device or a new one will then be returned to you. In the case of defects which occur after the end of the warranty period, the devices can also be returned. However, repairs which become necessary after the end of the warranty period will be subject to a service fee.

Important:

Make sure to return the device carefully packed in the original packaging in order to prevent transport damage. Please also enclose the cash receipt (or a copy). This warranty does not imply any restriction of your statutory rights.

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Un « Certificat de conformité » conforme aux directives applicables, établi par la société Bresser GmbH est disponible.